Panasonic



MDF-U33V-PA

VIP[®] Series

-86°C Ultra-Low Temperature Upright Freezer 11.8 cu.ft. 333 L



Panasonic's VIP Series ultra-low temperature freezers are designed for safe, reliable storage of biologicals. Energy efficiency in VIP Series freezers is achieved through balanced temperature performance and power management. The advanced combination of cabinet design, electronics, refrigeration and critical components are engineered for enhanced security, performance and reliability.

Reliability You Can Count On

The robust refrigeration system is powered by Panasonic engineered compressors. The system provides dependable cooling by balancing temperature performance and energy management. Reserve refrigeration power assures exceptional temperature recovery following door openings. VIP freezers have demonstrated >99% reliability with excellent performance history.

Maximum Storage, **Minimal Footprint**

The upright freezer employs a patented VIP insulation composite that maximizes space efficiency and holds up to 192 2" boxes or 144 3" boxes within a 6.25 ft² (0.58 m²) footprint.

Uniform **Preservation**

The combination of patented VIP vacuum insulation panels with Panasonic designed compressors, specific to ultra-low temperature applications, assures superior uniformity (+/- 5°C) throughout the cabinet to maintain the integrity of the samples.



Microprocessor Controls/Alarms

Integrated microprocessorbased controller and LED display include comprehensive setpoint, alarms, monitoring, diagnostic and communication features.*

*Control panel color & appearance may vary per equipment model.



Optimum Sample Storage

This space saving cabinet design offers a reduced wall thickness for maximum interior volume. Ideal for storage of DNA, RNA and other biological samples.



Extra Protection, Easy Access

The stainless steel framed, insulated inner doors with gasket seal firmly against the cabinet to help minimize change in interior temperatures during routine door openings. Inner door latches feature an ergonomic design for one-handed operation.



VIP -86°C ULT Upright Freezer




VIP Vacuum **Insulated Panels**

Inside-out engineering starts with a patented VIP insulation composite which results in an efficient, thin-wall cabinet. Insulated inner doors improve temperature uniformity, minimize cold air loss during door openings and extend warm-up times during power outages.

Model MDF-U33V-PA Temperature **Uniformity Data**


-80°C setpoint at 20°C ambient temperature, empty freezer

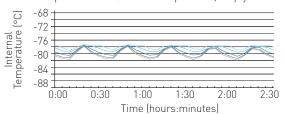

As interior air temperature stabilizes at the desired setpoint, mass storage volume sustains tight temperature uniformity of the load inside the insulated cabinet.





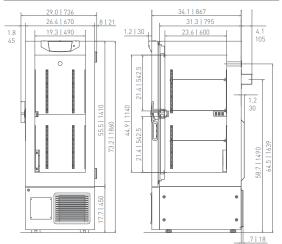

Panasonic

		MDF-U33V-F
Model Number		MDF-U33V-PA
External Dimensions (W × D × H) 11	inches mm	26.4 × 34.1 × 73.2 670 × 867 × 1860
Internal Dimensions (W \times D \times H)	inches mm	19.3 × 23.6 × 44.9 490 × 600 × 1140
Volume	cu.ft. liters	11.8 333
Net Weight	lbs. kg	562 255
Capacity (2" boxes)	qty	192
Capacity (3" boxes)	qty	144
Inventory Racks (2" boxes)	qty	12
Inventory Racks (3" boxes)	qty	12
Performance		
Warranty		Panasonic Five Year Warranty 2)
Cooling Performance 3	°C	-86
Temperature Setting Range	°C	-50 to -90 in 1°C increments
Temperature Control Range 3)	°C	-50 to -86 in 1°C increments
Control		
Controller		Microprocessor, touchpad data entry with display lockou
Display		LED
Temperature Sensor		Pt-1000
Refrigeration		
Refrigeration System		Cascade
Compressors	W	(1) 450 High-Stage and (1) 750 Low-Stage
Refrigerant	**	CFC-free R-404A and R-508A
Insulation Material		Rigid polyurethane foam (PUF) + VIP
Construction		Rigid polydrethane loant (FOF) + VIF
Exterior & Interior Material		Painted steel
Outer Door	qty	1
Outer Door Lock		Standard
Inner Doors	qty	2 insulated, ABS, w/ stainless steel frame and positive latches
Shelves	qty	3
Shelf Dimensions (W × D)	inches mm	18.3 × 21.1 464 × 535
Max. Load - per Shelf	lbs kg	110 50
Max. Load - Total Freezer	lbs kg	440 200
Vacuum Release Port		Standard
Access Ports	qty	2
Access Port Position		Center, upper back wall and back left chamber floor
Access Port Diameter	inches mm	0.6 17
Casters	qty	4 (2 leveling feet on front base)
Alarms		(V = Visual Alarm, B = Buzzer Alarm, R = Remote Alarm
Power Failure		V-B-R
High Temperature		V-B-R
Low Temperature		V-B-R
Filter		V-B
Door Open		V-B
Electrical and Noise Level		. 5
	A	115V, 1Ø, NEMA 5-20P,
Power Supply	•	requires NEMA 5-20R receptacle
Frequency	Hz	60
Noise Level ⁴⁾	dB (A)	49
Options		
Liquid CO₂ Back-up System		CKV-UB2I-PW
Temperature Recorder		
Circular Type	6" 7 day circular	MTR-C954 ⁵⁾
onar r upor	oz charto per box	57 1000001/LV
Ink Pen	pack of 6	R252
Circular Type Chart Paper	6", 7 day circular 52 charts per box	MTR-C954 ^{s)} C7100386REV

^{1]} Exterior dimensions of main cabinet only, excluding handle and other external projections

Wireless, cloud-based

LabAlert® Monitoring System

² Current warranty offered at time of printing and may be subject to change
³ Air temperature measured at freezer center, ambient temperature *30°C, no load
⁴ Nominal value - Background noise 20dB[A]
⁵ Requires sensor cover MTR-DU700SF-PW